



**Product Name:** HotShot™ Mastermix  
**Catalogue Number:** HS002  
Pack size: 5 x 1.0 ml  
Colour of lid: Blue  
Batch Number: 0306-1  
Expire Date: December 2008

**Double concentrated HotStart PCR mix, containing optimised mixture of *Taq* polymerase, anti-*Taq* polymerase monoclonal antibodies in 2 x Reaction Buffer (6 mM MgCl<sub>2</sub>) with 400 μM dNTPs and Stabiliser**

**For instant and accurate PCRs:**

- Thaw the **HotShot™ Mastermix**
- Take out the desired volume into the PCR tube.
- Add DNA, primers, fluorescence dye/probe and dH<sub>2</sub>O to an equal volume of the **HotShot™ Mastermix** (for instance to a 5ul volume of **HotShot™ Mastermix**, add 1ul DNA, 2ul Primers, 1ul dye and 1ul dH<sub>2</sub>O)
- Place in a Thermal Cycler
- Run cycling profile according to application

**Initial denaturation step at 95°C for 5 mins is needed to activate the *Taq* polymerase**

Refreeze the rest of the **HotShot™ Mastermix**. It can be frozen/thawed many times without loss of activity

Store at -20°C

For Research Only

P.O. Box 3059, Stourbridge, DY7 6YN, Tel: +44 845 644  
2368, Fax: +44 845 644 2367  
E-mail: [info@cadama.co.uk](mailto:info@cadama.co.uk)  
VAT registration No.: 803412082  
Company registration No.: 4561660



# Cadama Medical Ltd

Broaden Your Horizons

P.O. Box 3059, Stourbridge, DY7 6YN, Tel: +44 845 644  
2368, Fax: +44 845 644 2367  
E-mail: [info@cadama.co.uk](mailto:info@cadama.co.uk)  
VAT registration No.: 803412082  
Company registration No.: 4561660